

# Orion and Saturn packages

## Installation Instructions

### 1. Orion Package

#### 1.1. Full Package

Download the current version of the Orion package, whose name is in the form "OR\_xxx.rar" (please refer to the main WEB site for download via [www.OrionMicroscopy.com](http://www.OrionMicroscopy.com) or [www.SemImaging.com](http://www.SemImaging.com)).

Unzip the .rar archive in a temporary folder and run setup.exe. This will create a new folder containing the software and will create a new icon. Running the software without any Orion board in your computer will put the software in demo mode: all functions will be available except saving to your drive.

#### 1.2. Update files only

This zipped file contains only the critical files required to update your system. They are:

- The executable itself ("Orion32.exe") has to be placed in the folder of your current software version (refer to the icon properties to find which folder is involved);
- The two DLLs that have to be copied into your c:\Windows\System32 folder, overriding any other version if necessary.
- 

### 2. Saturn Package

Same procedure as above except that the name of the executable is "Saturn.exe" instead of Orion32.exe.